



Quick-Start Instructions

- Verify ventilation clearance:
 - REF12i/20i/25i: 10" (25.4 cm) above
 - REF45i, FZR20i/25i: 16" (40.6 cm) above
- Verify electrical – 120 V outlet
 - REF12i/20i/25i: 15A dedicated circuit
 - REF45i, FZR20i/25i: 20A dedicated circuit
- Install casters. Lock all front casters (see Manual)
- Adjust door tension and sag (see Adjust self-closing door in manual)
- Set up product simulation bottle(s) – 50% glycerine/50% water (refrigerator) or 60% glycerine/40% water (freezer) solution or equivalent
- REF/FZR20i/25i, REF45i – Open façade door to the right of the Touchscreen and connect the 9V battery to the chart recorder, if equipped
- Press power switch to turn unit ON
- The HIGH/LOW temperature alarms are pre-set to conform with NSF/ANSI 456 Standard for Vaccine Storage:
 - REF12i/20i/25i/45i: 46.4 F (8 C) / 35.6 F (2 C)
 - FZR20i/25i: 5 F (-15 C) / -50 F (-46 C)
- To alter alarm values, touch the Alarm icon () in the lower left corner of the screen. Change the high alarm by touching the number displayed in the box to the right of the HIGH label. Type your desired value in the keypad displayed to the right of the display box or use the up and down arrows to scroll to the desired temperature; press to store new value. Change the low alarm by touching the number displayed in the box to the right of the LOW label. Type your desired value in the keypad displayed to the right of the display box or use the UP and DOWN arrows to scroll to the desired temperature; press to store new value.
- If this is a keypad:
 - Default access code is 1-2-3-4-5-6 (Enter) (See operation manual for instructions to change access code)
- If desired change temperature display from Celsius to Fahrenheit
 - Press the Settings Icon (). Touch the box next to C or F.