

HEALTHCARE

The next level of control

The ultimate in overall reliability and high-performance temperature control, Follett's innovative Performance Plus undercounter refrigerators and freezers elevate critical product storage to a whole new level.

Engineered by the industry's proven leader in medical-grade products, the Performance Plus touchscreen control package puts every essential feature at your fingertips. From data collection and alarms to an electronic locking system, it's a snap to operate and is backed by Follett's long-standing commitment to superior service and support.



Touchscreen control package

Designed exclusively to enhance data management and product protection, the Performance Plus touchscreen control package offers value-added, next-generation features that are both smart and beautiful:

- Intuitive 7" capacitive touchscreen user interface
- Lighted cabinet
- Power loss/door open alarming functionality
- Lockable refrigeration settings
- Battery backup on alarm functions and data logging
- Data and event recording and on-screen display/graphing
- USB data export capabilities
- Optional touchscreen-integrated keypad access control


A home screen that takes you everywhere

Product temperature


Touch once to display the high/low temperature with time and date stamp

Touch twice to quickly reset the high and low temperatures

Touch three times to show the alarm set point display

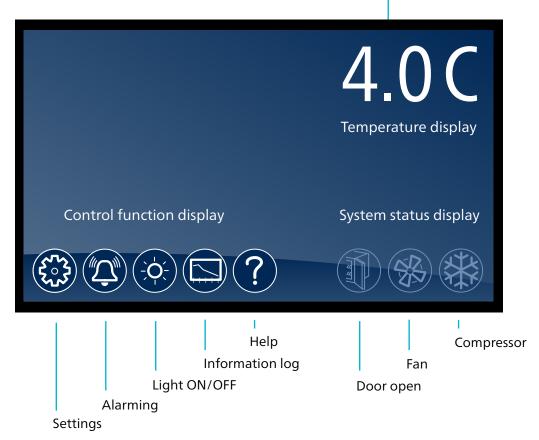

Easy access and complete control with just a touch

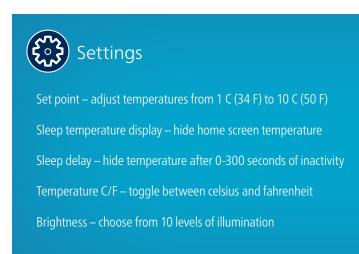





High and/or low alarm set points – select an alarm value

Audible alarm – toggle to engage/disengage audible alarms

Mute touch – mute audible alarm for up to 10 minutes





Easily access graphs, data and event logs

Zoom in or out on temperature graph

View all stored temperature data by simply scrolling



Data logging with just one touch

No need to drill down to access critical event and temperature information. Performance Plus puts everything you need just a screen touch away!


Data logging screen – easily access event and temperature logs



Event log – display a chronological listing of events, including alerts, alarms and errors



Temperature log – access a chronological listing of logged temperatures, starting with the most recent

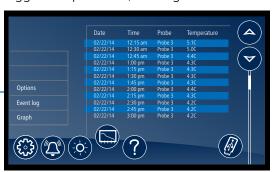

USB download



Saving event or temperature data has never been so easy!

- Simply insert a flash drive into the external USB port
- After the device is detected and confirmed, touch the ✓ in the yellow alert box to begin the data transfer directly from the event or temperature log screen
- Remove the device. It's just that simple.

Contextual help

Need help? Just touch the help icon located consistently at the bottom of each screen to access user-friendly answers and information.



System information

Touch the system info icon to display the model number, unit serial number and factory contact information without the hassle of searching for this information.



Optional onscreen keypad

Secure and straightforward, our optional keypad allows you to limit user access to your Performance Plus unit.

- Set code for authorized user entry
- Set up to 40 user codes
- Secure and reliable motor-driven bolt operation with spring-loaded latch
- Battery backup to safeguard the lock against power outages







FOLLETT HEALTHCARE is a trademark of Follett LLC.
FOLLETT is a registered trademark of Follett LLC, registered in the U.S.
Follett reserves the right to change specifications at any time without obligation.

